




High-Resolution Behavioral and Metabolic Phenotyping Systems for Rodents

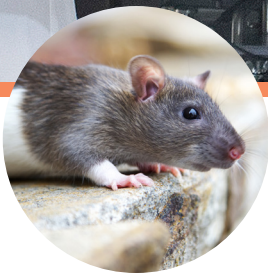

See What Others Miss



Promethion Core® is the only metabolic and behavioral phenotyping system that enables you to tightly synchronize metabolic data with behavioral events.

You will see every aspect of metabolism and behavior more clearly, with finer detail, quality, and repeatability—giving you complete confidence in your data.

Real-time data acquisition and analysis enables a new level of power to access, visualize, re-analyze and report study information. Sable's unique raw data storage protocol means there are no hidden algorithms or pre-conditioning of data—just fully traceable results that are available anytime for re-analysis. We make it easy for you to extract the parameters important to your research.

Promethion Core is designed by scientists, for scientists.


Our knowledgeable staff assists you every step of the way, including experimental design, system configuration, setup, training, and ongoing support.



Temporal Resolution Comparison: Promethion Core vs. Other Systems




Fast Response for Synchronized Results

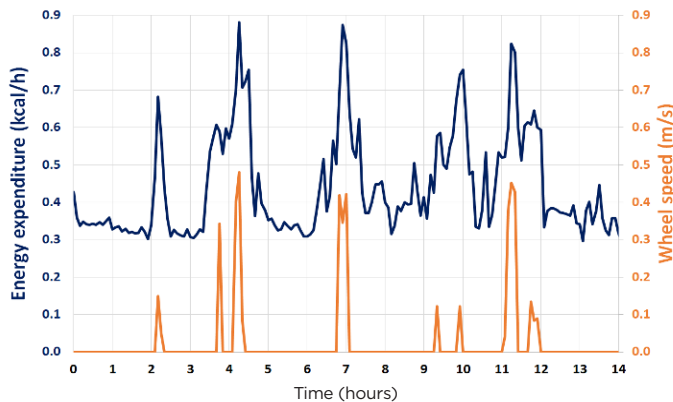

Promethion Core multiplexed system allows you to realize the full benefit of high resolution performance. The data above shows the synchronization of activity with energy expenditure. The EE graph shows sharp peaks and stable baselines enabling accurate active and resting energy expenditure calculation.

Competitor's System Slow, Blurred Response

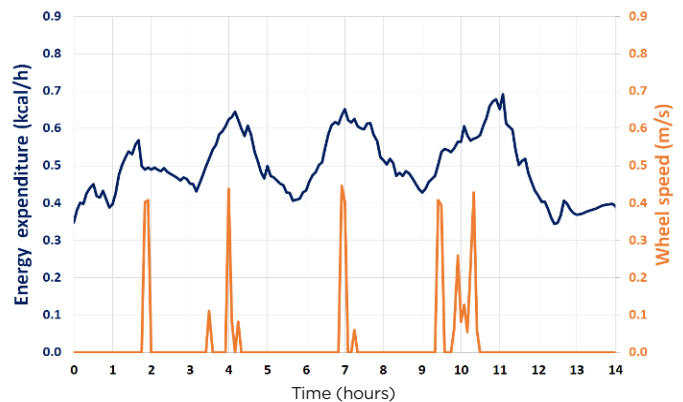

Competitors' continuous systems limit your research capabilities by using old, low resolution technology. The graph above shows broad peaks and no baseline. Therefore, the energy expenditure values are inaccurate.

Promethion Metabolic Treadmill

Promethion Core easily switches between metabolic cages and treadmills for controlled cardiometabolic research.

Fitness Testing: Exercise stimuli includes manual push broom and mouse-activated light, shock and auditory features.

Respirometry: Promethion Core system synchronizes changes in mouse metabolism with changes in treadmill speed and incline.

Automated Control: Hands-free incline and speed adjustments and automated self-leveling ensure incline accuracy and data fidelity.




Real-Time Monitoring: Promethion Live monitors and displays in real-time the respirometry measurements and the changes in incline and speed.

Animal Safety: The emergency stop button enables immediate access to the mouse. The shock grid is programmable not to exceed 5mA.

Real-Time Monitoring & Analysis

Setup with Confidence

The Promethion Live™ software platform maximizes experiment success by verifying system performance and simplifying experiment setup. One-step gas calibration ensures accurate respirometry measurements. User selectable configuration options are system specific. Status indicators of vital components verifies system performance. Routine experiment configurations are simply saved and reused. Smart setup delivers reliable experiments so that you can proceed with confidence.


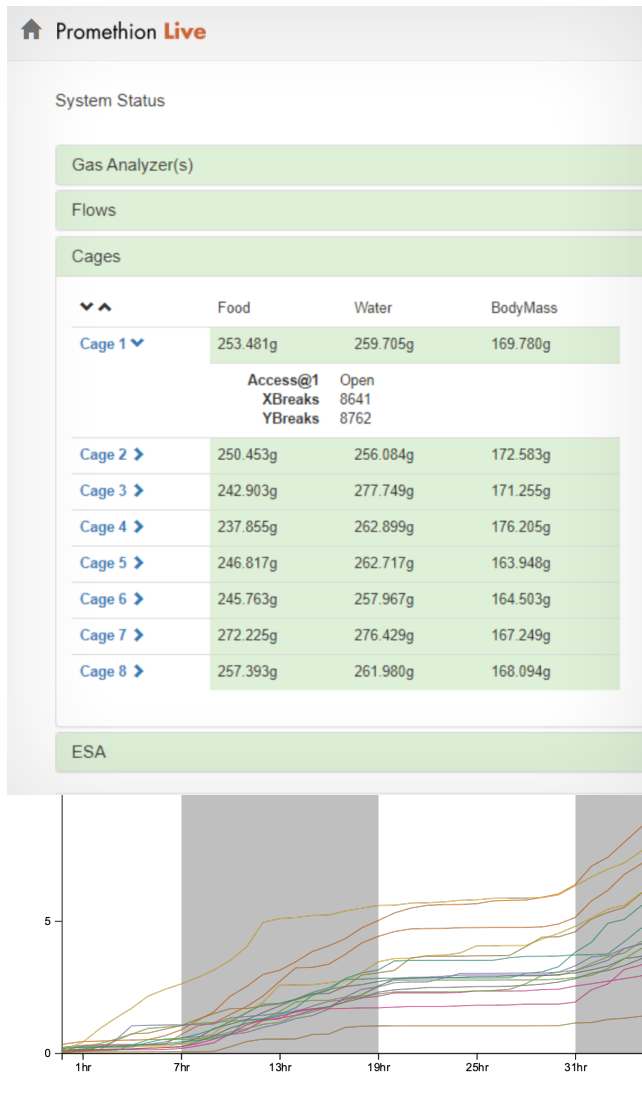


Actionable Intelligence

Promethion Live empowers you with real-time animal and system information needed to make key decisions at any time and any place. You'll always know the health and welfare of each animal. By monitoring activity, ingestive behavior and weight change, you will know when to make adjustments. You have total control over your experiment through continual monitoring of all system components.

Real Time Experiment Evaluation

Now you can explore comparative trends during an experiment looking at every facet measured, such as running activity, food consumption, or energy expended (data shown). The experiment result is no longer a black box to be opened at the end. Promethion Live provides the mean or median of grouped data for statistical comparison in real time. The data may be presented in real time or extracted for further evaluation.

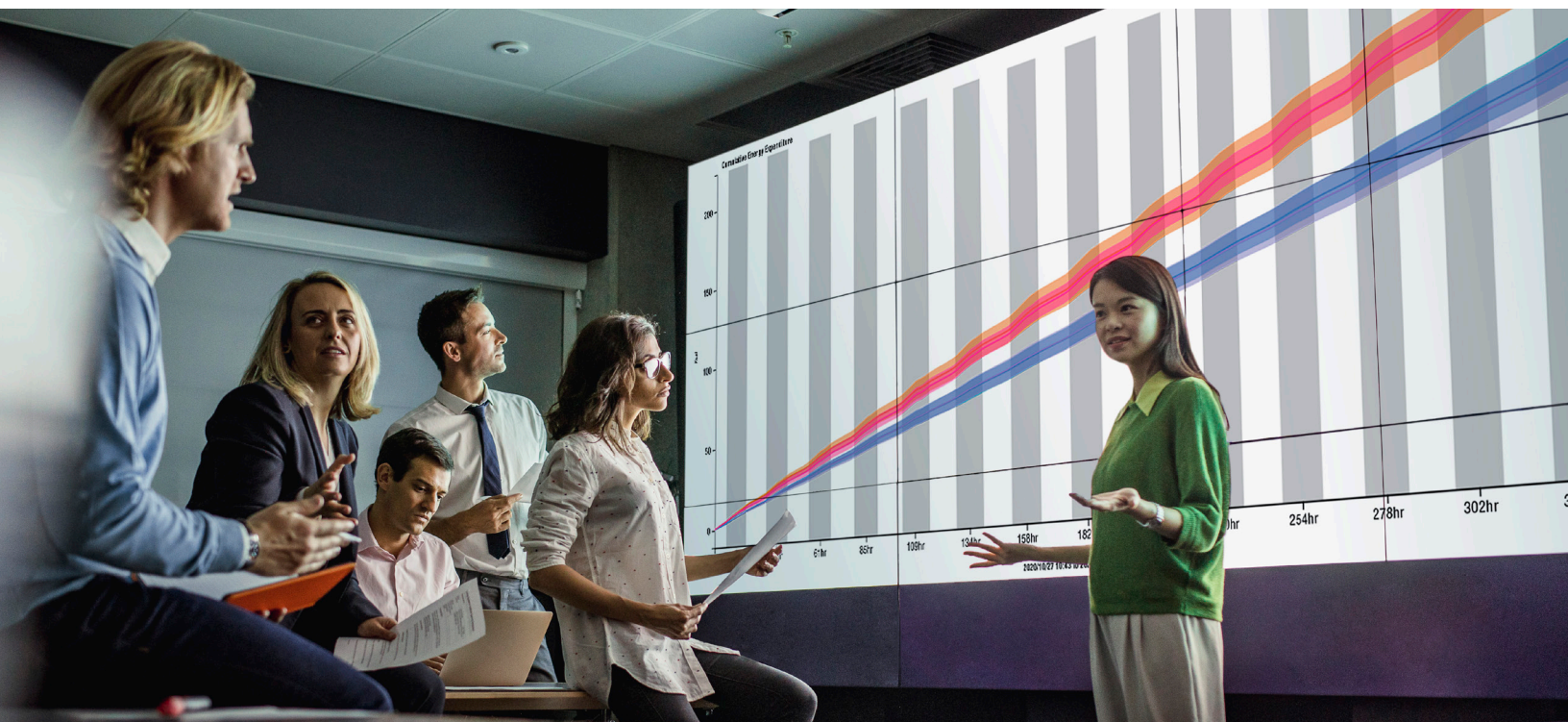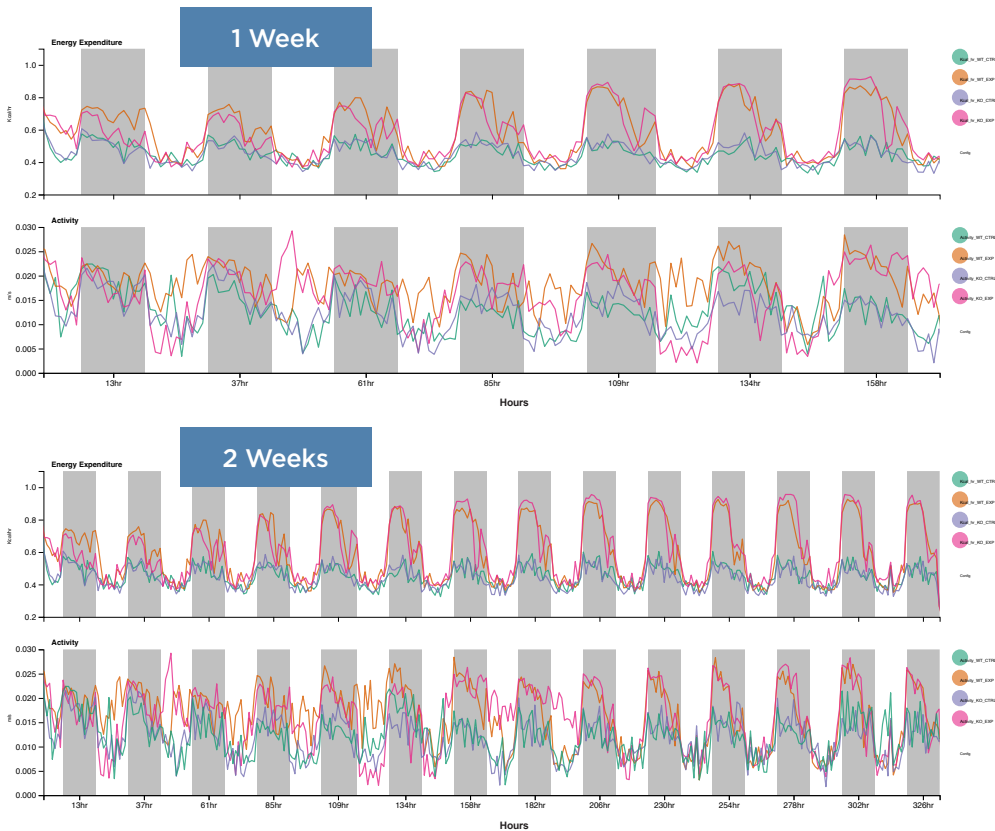

Metabolic Phenotyping Cage

Promethion Core systems are available in 8-16 cage configurations, with the ability to connect multiple systems for larger studies.

Exercise Research


The **Running Wheel Module** measures mouse running distance and speed. You can correlate running with metabolic measurements, including accurate active energy expenditure, and integrate running behavior into total behavior assessment.

Weight Assessment

The **Body Weight Monitor** measures weight changes in response to experiment or time.


Activity Monitoring

Promethion **Beam Break Activity Monitor** enables second by second analysis of total activity. This monitor can easily be used stand alone, or synchronized with metabolic measurement and other elements of the Promethion system.

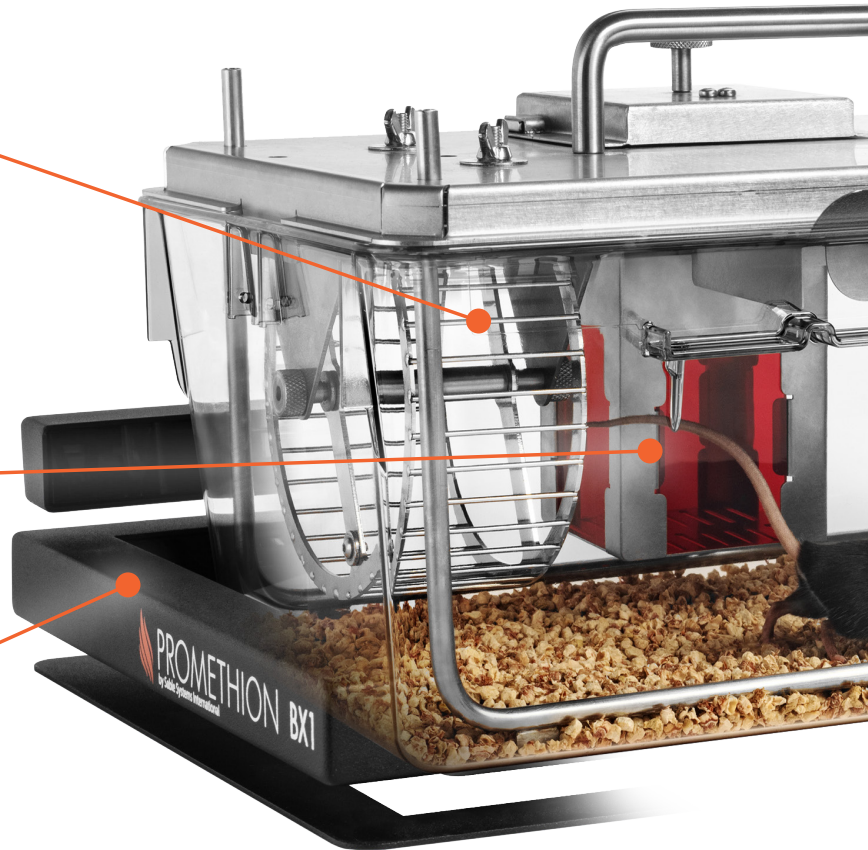


Gnotobiotic or Microbiome Research


The **Bioexclusion and Biocontainment Mouse Cage** is used for gnotobiotic or microbiome studies and includes 2-way HEPA filtration.

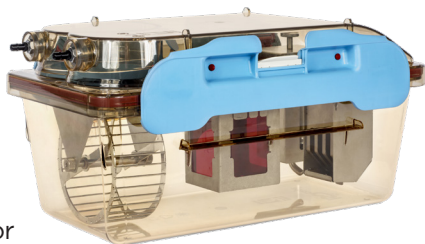

Rat Research

The **Rat Cage** has all the features and performance benefits of the standard mouse metabolic cage.

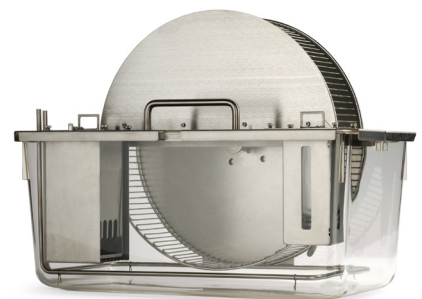

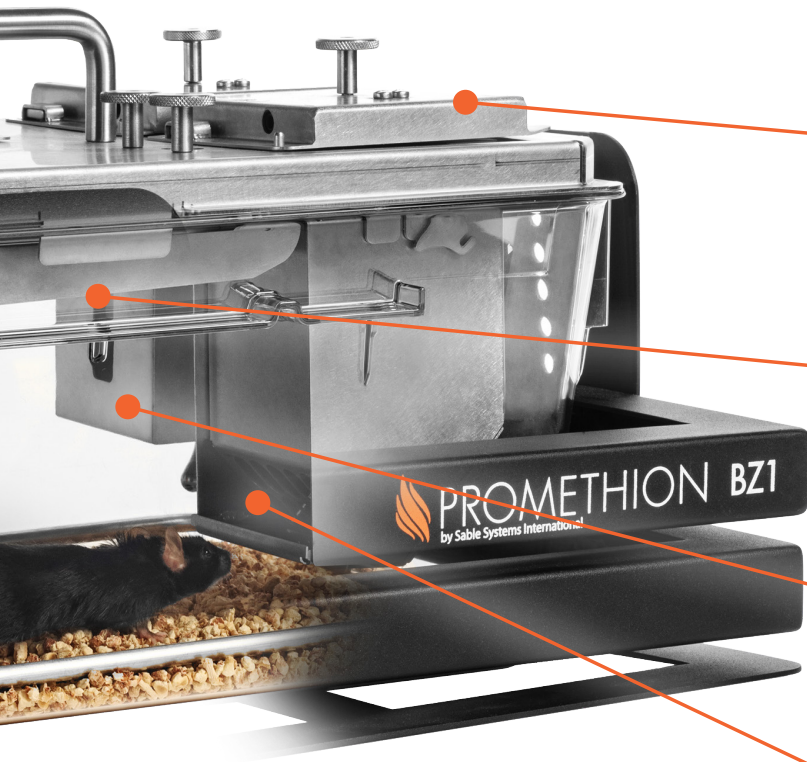

Fully Configured
Respirometry
Cage for Mice

Food and Water Handling

Magnetic Mounts simplify cage maintenance and workflow. They allow for quick swapping of the food, water and body mass measurement modules.

Ingestive Behavior Studies

The **Access Control Door** added to the food module enables caloric restriction, paired, and yoked feeding studies.

Water Intake Assessment


The **Water Intake Module** measures real-time liquid intake and allows you to conduct preference studies.

Food Intake Assessment

The **Food Intake Module** measures real-time food intake and allows you to conduct preference studies.


Ingestive Behavior
Research

The **Dual Access Mouse Cage** is designed for liquid or food preference studies and food access control studies.

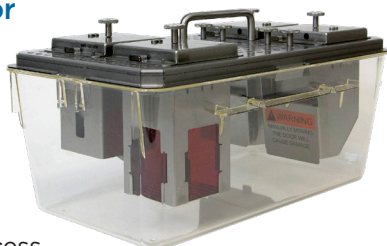

Rat Ingestive
Behavior Research

The **Group-House Rat Cage** measures intake events of individual rats. Designed for studies of liquid or food preference, and food access control.

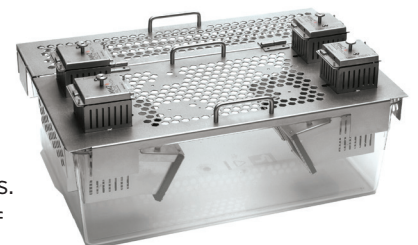

SABLE ENABLES

Promethion Core CGF

Promethion Core's Combined Gas & Flow (CGF) unit is a single organized housing that contains modular hot-swappable gas analyzers, flow regulators, pumps, onboard computer, and the Promethion Live software platform. This single integrated unit minimizes footprint and reduces clutter from tubing and cords.



Promethion Core Options


Environmental Control Cabinet

Perform temperature or light experiments on up to 16 animals in a single environmental cabinet specifically designed for rodent phenotyping studies.




Telemetry

Extend your research. Promethion Core couples with telemetry systems. Only Promethion Core has the resolution that enables correlation of metabolism data with the rapid signals produced by continuous glucose monitoring, pictured.

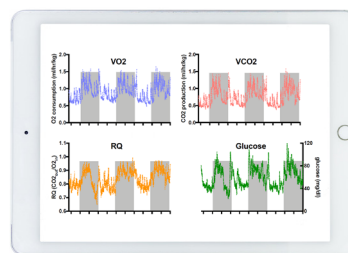


Stable Isotope Gas Analyzer

Study specific nutrient utilization by measuring ^{13}C or ^{18}O isotopes in exhaled breath.

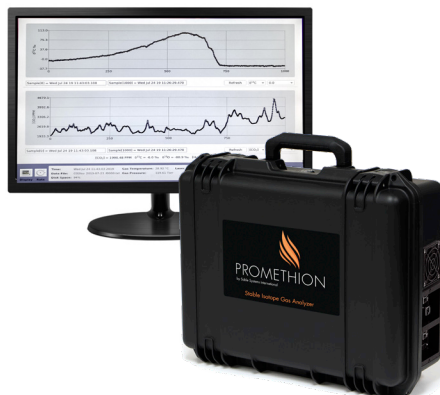

Methane Gas Analyzer

Study changes in gut, methanogenic microbes by measuring methane in exhaled breath.

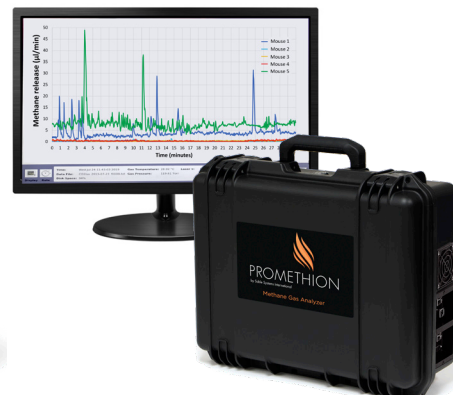

By Scientists, for Scientists



Sable Systems International
3840 N. Commerce Street
North Las Vegas, NV 89032, USA
TELEPHONE:
US: +1 800 330 0465 / +1 702 269 4445
EMAIL: sales@sablesys.com



Sable Systems Europe GmbH
Ostendstr. 25
D-12459 Berlin, Germany
TELEPHONE: +49 30 5304 1002
FAX: +49 30 5304 1003
EMAIL: sales@sablesys.eu

www.sablesys.com